4T1-Fluc-Neo/iRFP-Puro



Product Description

Product Name: 4T1-Fluc-Neo/iRFP-Puro

Catalog Number: CL078 Lot Number: CL-IM168

Species: Mouse (Mus musculus)

Strain: BALB/cfC3H

Cell type: Mammary carcinoma
Parental cells: 4T1 (ATCC® CRL-2539TM)*

Morphology: Epithelial Growth mode: Adherent

Reporter genes: Firefly luciferase (Fluc)

Near infrared fluorescent protein (iRFP)

Selection genes: Neomycin (Neo)

Puromycin (Puro)

This is a polyclonal population derived from the mammary carcinoma 4T1 cell line (ATCC® CRL-2539™). Parental 4T1 cells were transduced with 1) LV-Fluc-P2A-Neo (Imanis #LV011) encoding the firefly luciferase (Fluc) cDNA under the spleen focusforming virus (SFFV) promoter and linked to the neomycin resistance gene (Neo) via a P2A cleavage peptide and 2) LV-iRFP-P2A-Puro (Imanis #LV032) encoding the near infrared fluorescent protein (iRFP; ex./em = 690/713) cDNA under the SFFV promoter and linked to the puromycin resistance gene (Puro) via a P2A cleavage peptide. High Fluc and iRFP expressing cells were selected using G418 and puromycin. The lentiviral vectors are selfinactivating (SIN) vectors in which the viral enhancer and promoter have been deleted. Transcription inactivation of the LTR in the SIN provirus increases biosafety by preventing mobilization by replication competent viruses and enables regulated expression of the genes from the internal promoters without cis-acting effects of the LTR1.

*The ATCC trademark and any and all ATCC catalog numbers are trademarks of the American Type Culture Collection.

Mycoplasma Testing

This cell line has been tested for mycoplasma contamination and is certified mycoplasma free.

Cell Line Authentication

The parental 4T1 cell line was authenticated and certified free of interspecies cross contamination by STR profiling with 27 STR loci.

Recommended Uses

These cells are suitable for *in vitro* and *in vivo* experimentation.

The Fluc and iRFP transgenes facilitate *in vivo* noninvasive imaging of implanted cells.

Fluc and iRFP are immunogenic and may cause tumor rejection in immunocompetent mice. For the most consistent results, immunocompromised mice are recommended for studies.

References

¹Miyoshi et al. J Virol. 1998. 72:8150-8157.

Storage Instructions

Remove cells from the dry ice packaging and immediately store in the vapor phase above liquid nitrogen (below -130°C).

Complete Growth Medium

RPMI-1640 Medium (RPMI) 10% fetal bovine serum (FBS) 1% Penicillin/Streptomycin 0.1 mg/mL G418 2 µg/mL puromycin

G418 and puromycin should \underline{NOT} be added to the medium until a culture has been well established from the thawed cells (about 1 week). It is also recommended that a backup frozen cell stock be generated (see below) before adding G418 and puromycin to the growth medium.

Caution! Typical commercial puromycin stocks are provided at a concentration of 10 mg/mL or 10,000X.

Thawing Instructions

- 1. Thaw cells by gently swirling in a 37°C water bath. To limit contamination, do not submerge the O-ring and cap.
- 2. When cells are ~70% thawed (less than 1 min), remove the vial and wipe down with 70% ethanol. Allow tube to dry completely.
- 3. In a biosafety cabinet, transfer the cells into a 15 mL conical tube containing 5 mL of complete growth medium. Centrifuge cells at ~250 x q for 3-5 min.
- Remove supernatant and resuspend cells in 1 mL complete growth medium. Transfer cells to a T75 flask containing 10 mL complete growth medium.
- 5. Incubate the culture at 37°C with 5% CO₂. Cells should reach full confluency 2-3 days after thawing.

Subculturing Instructions

Volumes are given for a T75 flask. Increase or decrease as needed.

- 1. Remove culture medium from cells.
- Carefully wash the cell monolayer with 5-10 mL of phosphate buffered saline.
- Add 2 mL of 0.25% Trypsin-EDTA solution to the flask and incubate at room temperature until cells have dissociated (approx. 2-5 min).
- 4. Neutralize the trypsin by adding 8 mL complete growth medium and mix by gently pipetting up and down.
- Transfer desired portion of the cells to a fresh T75 flask. Add fresh complete growth medium to a total volume of 10 mL and return cells to 37°C/5% CO₂ incubator.

For maintenance, a subcultivation ratio of 1:10 is recommended. At this ratio cells will be ready for passage every 3-4 days. 4T1 cells frequently clump and should be passaged when they reach 80-90% confluency overall. Cell clumping is especially common after initial thawing and cells may need to be passaged more frequently at a lower subcultivation ratio for 2-3 passages after thawing.

Freezing Medium

Cells can be amplified and used to generate additional frozen stocks. Frozen stocks should be preserved in a designated cryopreservation medium or in complete growth medium without G418 and puromycin supplemented with 5-10% DMSO.

4T1-Fluc-Neo/iRFP-Puro




Certificate of Analysis

Testing performed by Imanis Life Sciences

Test description	Result
Post thaw viable cell recovery	98% viability
Sterility	No contamination detected
Mycoplasma	No contamination detected
Neomycin selection	Pass QC
Puromycin selection	Pass QC
Luciferase expression	Pass QC
Fluorescence expression	Pass QC
Average doubling time	14.6 h*


^{*}Doubling time represents the average doubling time during logarithmic growth. This value should be used for general estimation only.

Morphology:

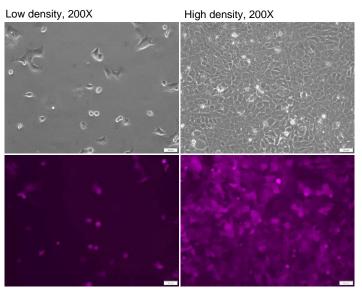


Low and high density photos taken after thawing.

Luciferase Expression:

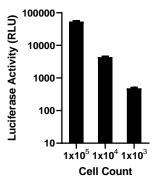

The indicated number of cells were placed in wells of a 96-well plate. After the addition of 15 mg/mL d-luciferin, bioluminescence was immediately read using a microplate reader.

Fluorescence Expression:

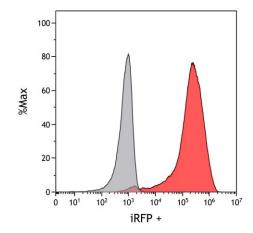

4T1-Fluc-Neo/iRFP-Puro (red) or isotype control (4T1; grey) cells were fixed with paraformaldehyde and analyzed by flow cytometry.

Legal Disclaimers

LIMITED PRODUCT WARRANTY
THIS WARRANTY LIMITS OUR LIABILITY TO REPLACEMENT OF THIS PRODUCT. NO OTHER WARRANTIES OF ANY KIND,
EXPRESS OR IMPLIED, INCLUDING, WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR
A PARTICULAR PURPOSE, ARE PROVIDED BY IMANIS. IMANIS SHALL HAVE NO LIABILITY FOR ANY DIRECT, INDIRECT,
CONSEQUENTIAL, OR INCIDENTAL DAMAGES ARISING OUT OF THE USE, THE RESULTS OF USE, OR THE INABILITY TO USE
THIS EDOPALIES.

FOR IN VITRO USE ONLY. THIS CERTIFICATE IS A DECLARATION OF ANALYSIS AT THE TIME OF MANUFACTURE.

PURCHASER NOTIFICATION

LIMITED LICENSE NOTICE - RESEARCH USE ONLY

THE IMANIS CELL LINES ARE NOT INTENDED FOR USE IN HUMANS. CELL LINES TRANSDUCED WITH LENTIVIRAL VECTORS ARE CLASSIFIED AS BIOSAFETY LEVEL 2 REAGENTS AND SHOULD BE USED UNDER THE APPROPRIATE BIOSAFETY

THE IMANIS CELL LINES ARE NOT INTENDED FOR USE IN HUMANS. CELL LINES TRANSDUCED WITH LENTIVIRAL VECTORS ARE CLASSIFIED AS BIOSAFETY LEVEL 2 REAGENTS AND SHOULD BE USED UNDER THE APPORIATE BIOSAFETY LEVEL PER INSTITUTIONAL GUIDELINES.

THE PURCHASER AGREES THAT IMANIS MATERIALS DESIGNATED AS BIO-SAFETY LEVEL 2 CONSTITUTE KNOWN PATHOGENS AND THAT OTHER IMANIS MATERIALS NOT SO DESIGNATED AND ANY PROGENY OR MODIFICATION MAY BE PATHOGENIC UNDER CERTAIN CONDITIONS. PURCHASER ASSUMES ALL RISK AND RESPONSIBILITY IN CONNECTION WITH THE RECEIPT, HANDLING, STORAGE, DISPOSAL, TRANSFER AND USE OF THE IMANIS MATERIALS INCLUDING WITHOUT LIMITATION TAKING ALL APPROPRIATE SAFETY AND HANDLING PRECAUTIONS TO MINIMIZE HEALT OR ENVIRONMENTAL RISK, PURCHASER AGREES THAT ANY ACTIVITY UNDERTAKEN WITH THE IMANIS MATERIALS AND ANY PROGENY OR MODIFICATION WILL BE CONDUCTED IN COMPLIANCE WITH ALL APPLICABLE QUIDELINES, LAWS AND REGULATIONS. THE IMANIS MATERIAL, ANY OTHER IMANIS PRODUCTS, AND ANY TECHNICAL INFORMATION AND ASSISTANCE PROVIDED BY MANIS ARE PROVIDED AS IS, WITHOUT WARRANTIES OF ANY KIND, EXPRESS OR IMPLIED, INCLIDING BUT NOT LIMITED TO ANY IMPLIED WARRANTIES OF MERICHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, MANUFACTURE ACCORDING TO GAMP STANDARDS, TYPICALITY, SAFETY, ACCURACY AND NON-INFRINGEMENT, IN O EVENT SHALL IMANIS, ITS PARENTS, SUBSIDIARIES, DIRECTORS, OFFICERS, AGENTS, EMPLOYEES, ASSIGNS, SUCCESSORS AND AFFILLATE (COLLECTIVELY "IMANIS INDEMNIFIED PARTIES") BE LIABLE FOR INDIRECT, SPECIAL, INCIDENTAL OR CONTRACT, TORT, NEGLIGENCE, STRICT LIABILITY, STATUFE OR OTHERWISS EVEN IF IMANIS HAS BEEN ADVISED, KNEW OR SHOULD HAVE KNOWN OF THE POSSIBILITY OF SUCH DAMAGES, INCLUDING, BUT NOT LIMITED TO, LOST PROFITS, CONTRACTOR OF SUBSTITUTE PROPOLETO'S OR CLAIMS OF LICABLES CUSTOMERS FOR SUCH DAMAGES, INCLUDING, BUT NOT LIMITED TO, LOST PROFITS, AGREEMENT FOR THE TWELVEL (2) MONTH PERIOD PRECEDING THE ACTUAL AMOUNTS PAID BY PURCHASER UNDER THIS AGREEMENT FOR THE TWELVEL (2) MONTH PERIOD PRECEDEN OF THE DATE OF THE EVENT GYN

Quality control by: AWD Quality Assurance by: RLV Effective Date: 16-Nov-2022